

Genie III v3.18.11 Software Update Install Instructions

1. Turn on and connect your Genie instrument to the computer.
2. Extract the .zip file to your computer and copy the files "GenieIII_Kernel_3_18_11_08Dec21.mot", "GenieIII_app_3_18_11_08Dec21.mot", and **if updating from versions prior to v3.18.2** "LoadSettings.set" and "UpdateWifi.set" and the **SYSTEM** folder to the Genie's drive **root directory**, replacing any files it asks about. This will take a few moments.
3. On the Genie, go to Settings, Utilities, and press Update and select "**GenieIII_Kernel_3_18_11_08Dec21.mot**" from the file list and press the tick button. The instrument will restart.
4. Go to Settings, Utilities and press Update and select "**GenieIII_app_3_18_11_08Dec21.mot**" from the file list and press the tick button. The instrument will restart.

If the software does not move off of the logo screen, hold down on the screen and a small pop-up menu will appear, select 'Install' then select "**GenieIII_app_3_18_11_08Dec21.mot**".

If it asks for a touchscreen calibration, do this and then try and hold down on the screen again. The instrument will again restart after installing the file.

5. Touch the screen again. If it asks "**Enter Manufacturer mode**". **Press No to this.**
6. **If updating from versions prior to v3.18.2**, go to Settings, Utilities, and press update. Select "**LoadSettings.set**".
7. Restart the instrument by using the soft power button in the bottom left corner of the screen.
8. **If the instrument is a MODEL: GEN3-02-02 ONLY and If updating from versions prior to v3.18.2** Go to Settings, Utilities, and press update. Select "**UpdateWifi.set**".
9. Restart the instrument by using the soft power button in the bottom left corner of the screen.

The instrument is now ready to use.